

A GUIDE TO GOWNING

1 GLOVES (ISO 3+) OR GLOVE LINERS (ISO 6+)



Select gloves of proper size and don them. Two layers of gloves ensures no contamination if top layer rips or tears.

2 FACEMASK (ISO 3+)



Select proper size and bend the nose-piece to ensure a snug facial fit. Check for leakage; nose and mouth should be covered

3 HOOD (ISO 3+) OR HAIR COVER (ISO 6+)



For hood, fully adjust to comfortable fit with the correct seal around the face area. For hair cover make sure whole scalp is covered.

4 GOGGLES (ISO 3+)



Wear goggles to cover the exposed eye area. Make sure goggles are placed firmly on the face and, if available, use adjusters on the strap to fasten accordingly.

5 SOCKS (ISO 3+)



Remove personal socks first and then remove cleanroom socks from packaging. Place over foot making sure sock is on securely

6 COVERALL (ISO 3+) OR LAB COAT (ISO 6+)




Step into the coverall or put on lab coat, ensure upper garment and sleeves do not touch the floor. Tuck hood fully inside the collar.

7 OVERBOOTS (ISO 3+) OR SHOE COVERS (ISO 6+)



Place over your shoes making sure they are fully covered from toe to heel. If wearing overboots, make sure they are fully fastened.

8 GLOVES (ISO 3+)

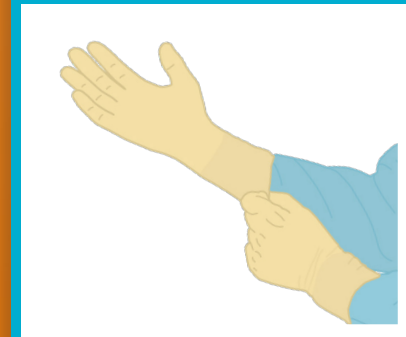

Follow the glove gowning procedure, don gloves and ensure the cuffs cover the coverall sleeve. You are now fully gowning for the cleanroom